Alicyclobacillus acidoterrestris- Guaiacol-Ident Kit

Reagents for the detection of *A. acidoterrestris* and potentially guaiacol producing *Alicyclobacillus* spp.
English: page 3 - 11

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***Attention:**

This kit contains only reagents used for amplification.
For the detection of amplicates the following additional kit is required:
Milenia® GenLine PCR-Universal module MGUP 1 (50 tests)

Table of Contents

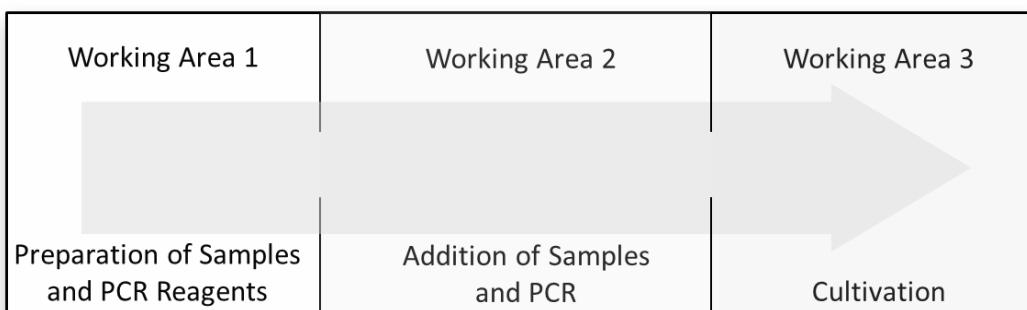
| | |
|---|----|
| Explanation of Symbols | 3 |
| Warnings and Precautions | 3 |
| Materials supplied, Storage and Stability | 4 |
| Materials Required (not included in the test kit) | 4 |
| Intended Use | 4 |
| Method and Test Principle | 4 |
| Procedure of the Test – A short Overview | 5 |
| Preparation of Samples and PCR Reagents | 6 |
| Test Procedure | 8 |
| Interpretation of Results | 9 |
| Trouble shooting | 10 |
| Literature | 10 |

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Warnings and Precautions

- Read this instruction manual carefully before running the test. Follow the sequence of the work steps
- Pay attention to the expiration dates of the individual components of the test
- Complete thawing of all components prior use is required
- Do not exchange components with different Lot numbers
- Store all reagents at **-15 °C to -20 °C** in the original packaging
- **Immediately after use**, store all components at **-15 °C to -20 °C**
- The disposal of waste materials must be carried out according to current local regulations
- Wear laboratory coat and protective gloves
- In order to avoid contaminations, run the different steps of the test at different locations
- The place where the reaction mixes are prepared must be separated from the place where cultures are handled.
- Decontamination of the work area and the equipment used must be carried out regularly!

Materials supplied, Storage and Stability

| REF: MGAaG Z | | | | | | |
|--------------------------------|----------|----------|---------|--------------|------------------|-------------------|
| Component | Part-Nr. | Quantity | Content | Preparation | Storage | Shelf life |
| PCR-Polymerase Master-Mix (2x) | MGPOL | 2 | 330 µL | Ready to use | -15 °C to -20 °C | until expiry date |
| PCR- Primer mix (2x) | MGPAaG | 2 | 280 µL | Ready to use | -15 °C to -20 °C | until expiry date |
| Positive Kit-Control | MGAaGPC | 1 | 50 µL | Ready to use | -15 °C to -20 °C | until expiry date |
| PCR-Water* | MGH2O | 1 | 1 mL | Ready to use | -15 °C to -20 °C | until expiry date |

*PCR-Water (MGH2O) can be used as diluent or negative control.

Material safety data sheets are available on request or from the company website
(www.milenia-biotec.de)

Materials Required (not included)

Important:

In addition to the components supplied with the kit, the following reagents and items are required:

- **Milenia® GenLine PCR-Universal module (MGUP 1).** Contains test strips and chase buffer.
- **Extraction Buffer (MGEXBA)**
- **Thermocycler Labcycler 48s**, Order code: MLCY, available at Milenia Biotec GmbH
- **Deep freezer:** -20 °C
- Pipets 2- 20 µL, 20- 200 µL, 100- 1,000 µL
- Sterile (DNA-free) pipet tips (with contamination protection for PCR)
- Sterile (DNA-free) reaction tubes 1.5 mL
- Sterile (DNA-free) PCR-tubes 0.2 mL e.g.: Supplier: Sorensen Bioscience Inc.
- 96-well microtiter plate
- (Desktop) centrifuge for reaction- and PCR-tubes
- Vortex mixer
- Magnetic stirrer
- Tube holder for reaction- and PCR-tubes, e.g.: isofreeze PCR tube rack

Intended Use

The Milenia® GenLine *Alicyclobacillus acidoterrestris-Guaiacol-Screen* is designed for the qualitative detection of *Alicyclobacillus acidoterrestris* and potentially guaiacol producing organism. This test follows if the *Alicyclobacillus* spp.-screen (**MGScACB 1, has to be ordered separately**) is positive.

Method and Test Principle

Milenia® GenLine tests are rapid detection systems that combine PCR (Polymerase Chain Reaction) methodology with detection and interpretation of the resulting products by lateral flow tests.

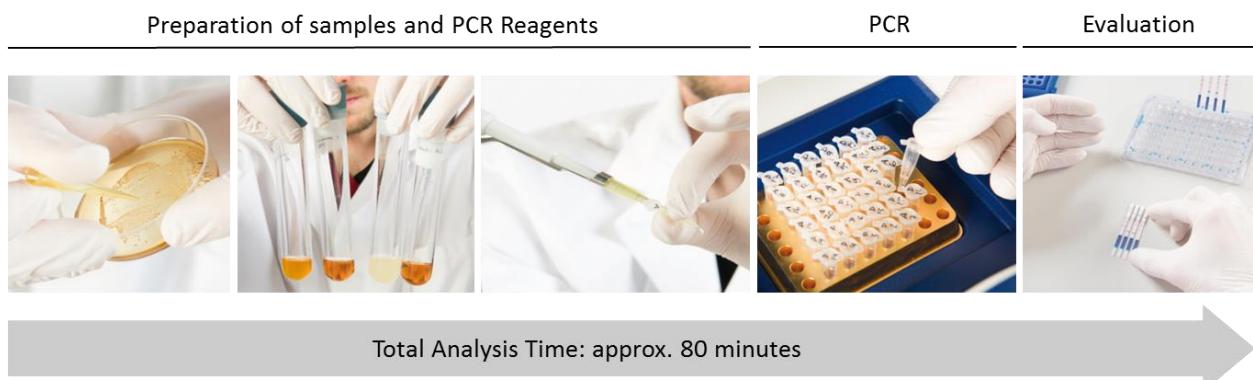
In a first part of the test, labelled specific primers (short, single stranded DNA) are used to amplify specific DNA fragments. In this process the labels are incorporated in the resulting PCR products and can be detected by a lateral flow test strip in a second step.

By the pre-performed *Alicyclobacillus* spp.-screen it was shown that the PCR process works properly. An additional amplificat control is not necessary.

Procedure of the Test – A short Overview

The entire procedure of the test can be divided into 3 different steps.

1. The PCR reagents and the samples are prepared.
2. After the addition of the sample to the PCR reagents, the PCR is started.
3. The resulting PCR products are detected by a simple lateral flow test strip.

Preparation of Samples and PCR Reagents

The preparation of samples and reagents includes the preparation of the PCR reaction mix and the samples to be analysed.

A) Sample Preparation

Samples to be analyzed are fruit juice concentrates, which were 1:10 diluted in BAT-B (Döhler GmbH, Darmstadt), heat activated and cultivated over a maximum of 5 days (according to IFU12) to enrich potentially existing *Alicyclobacillus*. DNA extract from sample preparation for *Alicyclobacillus* spp. screen should be already available. In this case use 2 µL out of the separately stored DNA-containing supernatant. Follow the instruction **if no DNA-extract** is available to setup the PCR.

Procedure:

- 1) Add 800 µl sterile water and 200 µl of the enriched probe in a 1.5 mL reaction tube, mix well and centrifuge for 3 min at 16 000 xg.
- 2) Homogenise the extraction buffer (MGEXBA) with the magnetic stir bar. Discard the supernatant from step 1) and add 200 µl extraction buffer (MGEXBA) to the pellet. Subsequently, the pellet is completely resuspended by thoroughly vortexing for minimum 15 s.
- 3) Incubate the suspension for 10 min at 95°C and mix thoroughly for 15 s (vortex).
- 4) Centrifuge the suspension for 1 min at 16.000 xg.
- 5) Carefully collect 2 µl of the supernatant and use in PCR. Please avoid the transfer of pieces from the pellet into the PCR reaction tube!

For later analysis the **pure** supernatant (without pellet) can be transferred into a new reaction tube and stored by -20°C for several months.

Analysis of single colonies from solid media (z.B.: BAT-A, Döhler GmbH):

Procedure:

- 1) Touch a colony on the agar plate with a pipet tip. Do not wipe across the colony.
- 2) Resuspend the cells in 100 µL sterile Water.
- 3) Add 2 µL of the cell suspension to the PCR reaction mix.

If too many cells are added to the mix, this may have a negative impact on the test. Therefore, a careful light touch in step 1) is recommended.

B) Preparation of PCR reaction mix

Each individual tube of reaction mix has a total volume of 23 µL. In order to prepare the mix, the PCR-Polymerase Master Mix (2x) and the PCR-Primer mix (2x) are mixed in a ratio as given in the table below. Finally the mix is filled in 23 µL aliquots into sterile PCR tubes.

The precise and contamination free preparation of the mix is the prerequisite of a proper function of the test. It is recommended to label all reaction tubes prior pipetting the samples. The final and aliquoted reaction mix must be used directly and placed cool (e.g.on ice) during the addition of samples. In case a positive (MGAaGPC) and a negative (MGH2O) Control is included in the run, add 2 additional tubes to the number of samples introduced in the test run.

| Number of PCR reaction tubes | PCR-Polymerase Master Mix (2x) (Volume to be pipetted) | PCR-Primer mix (2x) (Volume to be pipetted) |
|------------------------------|---|--|
| 1 | 13 µL | 11 µL |
| 5 | 65 µL | 55 µL |
| 10 | 130 µL | 110 µL |
| 20 | 260 µL | 220 µL |

Procedure:

- 1) According to the number of samples (plus a positive and negative control) PCR-Polymerase Master Mix (2x) is pipetted into a sterile 1.5 mL reaction tube (volume is given in the table).
- 2) PCR-Primer mix (2x) is added to the tube (see table).
- 3) The mix of both components is mixed well (pipet, Vortex).
- 4) Take an appropriate number of sterile 0.2 mL PCR-reaction tubes.
- 5) Pipet 23 µL of mix into each PCR reaction tube.
- 6) Close the PCR-reaction tubes immediately and add the sample at a different location.

Test Procedure

The test procedure describes the start and termination of the PCR reaction as well as the subsequent lateral flow based detection of PCR products by a universal test strip (Milenia® GenLine Universal Module MGUP1).

C) PCR reaction

Procedure:

- 1) After addition of the sample to the PCR tube containing the reaction mix, close the tube and mix well.
- 2) Place all PCR tubes into a minicentrifuge and centrifuge the samples briefly.
- 3) All liquid should be free of air bubbles and located on the bottom of the PCR tubes.
- 4) Place the PCR tubes into a thermocycler.
- 5) Close the lid of the thermocycler.
- 6) Start the specific ACB/AaG program.
- 7) After completion of the program, stop the program and remove the PCR tubes from the thermocycler.
- 8) When the program is finished, the thermocycler holds a temperature of 8°C. PCR products must be stored in a refugator (app. 4°C) and shoulb be analysed subsequently (max. 24 h later).

D) Detection of PCR products

For the detection of PCR products the **Milenia® GenLine PCR-Universal Modul MGUP 1** is required.

It is strongly recommended to use pipet tips with filter and separate pipets for the addition of the sample to be analyzed and the evalutation of PCR products to minimize the risk of DNA contamination. Regular decontamination of the pipets and a separation of the different working steps at different locations are recommended.

Procedure:

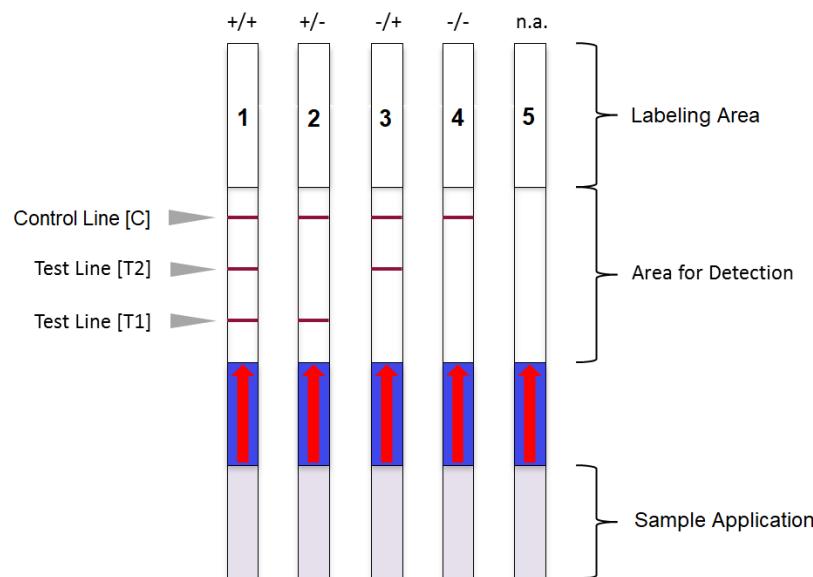
- 1) According to the number of tests, pipet 80 µL Chase buffer (MGCBA) each into clean wells of a 96 well microtiter plate.
- 2) Label each test strip (MGDS2A) needed for the analysis and take it aside.
- 3) Pipet 2 µL of each PCR product onto the sample application pad of a test strip.

- 4) Place the test strip with the sample application pad down into the Chase buffer of the wells of the microtiter plate and incubate for 5 min at room temperature. Read the results visually after exactly **5 min**.
- 5) For the documentation of the results, strips can be fixed onto a data sheet (MGDBES). An additional documentation by taking an image of the strips with a camera is recommended.

Results must be read exactly after 5 minutes! After prolonged time unspecific, weak signals may occur on the test strip which are not associated with positive results.

Interpretation of Results

Tests are interpreted visually. Potential test results are shown in the figure below.

The control line [C] confirms the proper function of the test strip. The test lines [T1] and [T2] are associated with the presence or absence of the specific target gene.

A) Positive Result (test strip 1, 2 and 3)

The prerequisite for a valid positive result is the appearance of a distinct test line [T1] and/or [T2], (test strip 1, 2 and 3).

The test line [T2] indicates whether it is *A. acidoterrestris*.

The test line [T1] indicates a potential guaiacol former.

The samples to be analyzed must first be tested with the *Alicyclobacillus spp.*-screen in order to exclude inhibitory effects in the PCR.

B) Negative Result (test strip 4)

A valid negative result requires a positive result in the *Alicyclobacillus spp.*-screen. This confirms that the PCR has worked properly. The test is negative, if the test lines [T1] and [T2] are not visible (Test Strip 4).

C) No analysis possible (test strip 5)

If no line is visible on the test strip, please contact the customer service at Milenia Biotec GmbH.

Trouble shooting

The negative control gives a positive result

Possible cause: DNA-Contaminations

Contaminations are a high risk when working with nucleic acids. Therefore the following preventive actions are strongly recommended:

- Check that the chosen PCR protocol of the thermocycler was correct.
- Decontaminate the workspace where the preparation of the reaction mix is done.
- Decontaminate the pipets.
- Use unopened, sterile plastic materials (Pipet tips/ 1.5 mL reaction vials/PCR vials)

After decontamination, a test run should be performed in order to check the effectiveness of the preventive actions.

It is recommended to run a positive and a negative control. The negative control **must** show a negative result.

If the negative control shows a positive result, this may be due to a contamination of the reagents. In this case new reagents (PCR-Water, PCR-Polymerase Master Mix (2x) and PCR-Primer Mix (2x)) must be used. In the subsequent control run, the negative control (PCR-Water, MGH2O) must be negative.

If the negative result shows a positive result again, please contact the customer service at Milenia Biotec GmbH.

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IFU: Method on the detection of taint producing *Alicyclobacillus* in fruit juices. IFU Handbook Microbiological Methods 3 (2004)

Additional Products available

| Product Name | Order No. | Content | Description |
|---|-----------|----------|--|
| Milenia® <i>Alicyclobacillus spp.</i> -Screen | MGScACB* | 48 tests | Reagents for the detection of <i>Alicyclobacillus spp.</i> |

For any questions and suggestions, please contact us:

Telefon: 0 641 – 94 88 83 – 0
Fax: 0 641 – 94 88 83 – 80
E-mail: info@milenia-biotec.de
Web: www.milenia-biotec.de

Alicyclobacillus acidoterrestris-Guaiacol-Ident Kit

Reagents for the detection of *A. acidoterrestris* and potentially guaiacol producing organism
English: Page 3 - 11

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Sind keine Linien auf dem Teststreifen zu erkennen, kontaktieren Sie bitte den Kundenservice der Milenia Biotec GmbH.
Ist die negative Kontrolle erneut positiv, kontaktieren Sie den Kundenservice der Milenia Biotec GmbH.