Mouse-Isotyping Sample buffers

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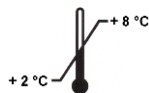
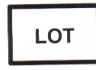
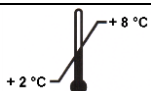
*** Attention:**

These reagents are designed for the application in Milenia[®] QuickLine Mouse Immunoglobulins-universal module (MQUM 1)

Table of Contents

Explanation of Symbols	3
Warnings and Precautions	3
Materials supplied, Storage and Stability	3
Materials Required (not included)	4
Intended Use	4
Principle of the Test	4
Test Procedure	4
Interpretation of Results	5
Trouble shooting	5
Assay Sensitivity	5
Additional available Products	5

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Warnings and Precautions

- Read this instruction manual carefully before performing the assay. Follow the sequence of the work steps
- Store reagents at **2°C to 8°C** in original packaging
- Check the expiration date printed on the vial and on each Sample buffer
- The disposal of waste materials must be carried out according to current local regulations
- Wear lab coat and protective gloves
- Material safety data sheets are available on request or can be downloaded from www.milenia-biotec.de

Materials supplied, Storage and Stability

REF: MQSBXX					
Component	Cat.-No.	Content	Preparation	Storage	Shelf life
Anti-mouse-IgG1 Sample buffer	MQSBG1	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse-IgG2a Sample buffer	MQSBG2A	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse-IgG2b Sample buffer	MQSBG2B	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse-IgG2c Sample buffer	MQSBG2C	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse-IgG3 Sample buffer	MQSBG3	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse-IgM Sample buffer	MQSBM	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse-IgA Sample buffer	MQSBA	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse-kappa Sample buffer	MQSBKLC	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse-lambda Sample buffer	MQSBLLC	0,5 mL	ready to use	2°C to 8 °C	until expiry date

Materials Required (not included)

Important:

- **Milenia[®] QuickLine Mouse Immunoglobulins-Universal module (MQUM 1)**
contains test strips for 100 tests
- **96-well microtiter plate**
- **Pipets (2-20 µL, 20-200 µL) and tips**

Intended Use

Milenia[®] QuickLine Mouse-Isotyping Sample buffers are designed for fast and easy identification of mouse antibody isotypes, subtypes and light chains in combination with the universal test strip (Milenia[®] QuickLine Mouse Immunoglobulins-Universal module (MQUM 1)).
For research use only. Not for diagnostic purposes.

Principle of the Test

Knowing the antibody-subclass is important when mouse monoclonal antibodies are produced. Hereby the optimal purification method can be chosen for the specific isotyp which results in the highest yield of the antibody. Furthermore positive clones can easily be detected with this test after recloning, if there is no antigen available.

Milenia[®] QuickLine Mouse-Isotyping is a lateral flow immunoassay designed for the determination of mouse monoclonal antibody isotypes, subtypes and light chains within 5 minutes.

The immunoassay can be used for cell culture supernatants and/or purified antibodies.

First the isotype-specific Sample buffer (e.g. anti-mouse IgG1) is pipetted into a well of a 96-well microtiter plate. Subsequently the test sample of unknown antibody-subclass is added. If a complex is formed between the Sample buffer and the test sample, binding to an anti-mouse antibody on the universal mouse test strip will result in a red line formation.

Test Procedure

Important notes:

All reagents must be equilibrated to room temperature (18-28°) before use.

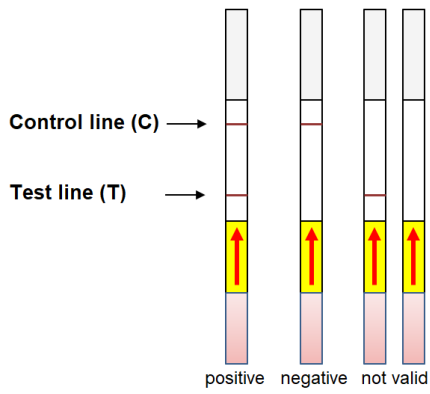
If the antibody concentration is too high, it can lead to a false negative result (high dose hook effect). This may occur when highly concentrated purified antibodies are applied in the test.

In this case, samples should be prediluted with 1 X PBS to enter in the working range of the assay (1 – 20 µg/mL).

Test setup:

1. Pipet 45 µL of the appropriate isotype-specific Sample buffer into the wells of a 96 well microtiter plate
2. Add 10 µL of the antibody test solution (about 1-20 µg/mL) and mix well
3. Place one universal dipstick of the **Milenia[®] QuickLine Mouse Immunoglobulins-Universal module** into every well
4. Incubate for **5 minutes** at room temperature
5. Read the results by visual inspection immediately after incubation time, test lines developed after this time are without relevance

Interpretation of Results

Positive: test line (T) and control line (C) are visible
Specific isotype has been **detected**

Note: faint red test line should be reported as positive

Negative: only control line (C) is visible
No isotype can be detected (valid negative result)

Not valid: only test line (T) or no line visible

Tests are interpreted visually. Examples of test results are shown in the figure above. The control line (C) should be positive in all cases, indicating a successful test. The test line (T) is associated with the presence or absence of the target isotype.

Trouble shooting

Problem	Possible cause(s)	Recommendation
No visible control line	Damaged dipstick	Repeat test with new dipstick
False negative result	a) concentration of purified antibody is too high b) concentration of antibody in cell culture supernatant is too low	a) dilute antibody in 1 X PBS b) use higher amount of antibody
Several test lines occur	a) Incubation time was exceeded b) Concentration of purified antibody is too high	a) Do not incubate for more than 5 Minutes b) Dilute antibody in 1 X PBS

If no line is visible on the dipstick please contact the Milenia Biotec GmbH customer service.

Assay Sensitivity

Sensitivity varies based on different conditions (purified antibody/cell culture supernatant). Antibody amounts between 1 and 20 µg/mL are detectable.

Additional Products available

Product Name	Order No.	Content	Description
Milenia® QuickLine Mouse Immunoglobulins-Universal module	MQUM 1	100 tests	Universal dipstick for mouse monoclonal antibody subtype detection

Mouse-Isotyping Sample buffers

Lateral flow reagents for the identification of mouse monoclonal antibody isotypes, subtypes and light chains
Only for use in combination with MQUM 1
English: Page 3-11

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Ist auf dem Dipstick keine Linie sichtbar, wenden Sie sich an den Kundendienst der Milenia Biotec GmbH.