milenia biotec

Mouse Antibody Isotyping Instruction

**Instruction for the Identification of Mouse Monoclonal
Antibody Iso-, Subtypes and Light Chains**

Required components:

Mouse Ab Isotyping Universal Module (Order No. MQUM 1)

Mouse Ab Isotyping Sample Buffers (Order No. MQSBXX)

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Web: www.milenia-biotec.de

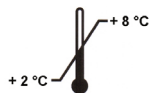
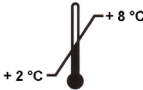
Table of Contents

Explanation of Symbols	3
Warnings and Precautions	3
Required Products, Storage and Stability	4
Additionally Required Materials	5
Intended Use	5
Principle of the Test	5
Test Procedure	5/6
Interpretation of Results	6
Troubleshooting	7
Assay Sensitivity	7
Documentation of Results	7
Quick Reference Guide	8

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Warnings and Precautions

- Read this instruction manual carefully before performing the assay. Follow the sequence of the work steps
- Store reagents at **2°C to 8°C** in original packaging
- Check the expiration date printed on the components
- The disposal of waste materials must be carried out according to current local regulations
- Wear lab coat and protective gloves
- Material safety data sheets are available on request or can be downloaded from www.milenia-biotec.de

Required Products, Storage and Stability

Isotyping monoclonal Mouse-Antibodies

Milenia QuickLine
Mouse Antibody Isotyping
Universal Module *1
(Order No. MQUM 1)

Lateral-Flow Universal Test Strips
essential for the test
2 x 50 tests

Milenia QuickLine
Mouse Antibody Isotyping
Sample Buffer *2
(Order No. MQSBXX)

9 Anti Isotype Buffers available
freely selectable and combinable
each Buffer 10 tests

*1 Milenia QuickLine Mouse Antibody Isotyping **Universal Module** (Order No. MQUM 1):

Product Name	Order No.	Content	Preparation	Storage	Shelf life
Universal Module	MQUM 1	2 x 50 test strips	ready to use	2°C to 8 °C Container must be always closed (protected from moisture)!	until expiry date

*2 Milenia QuickLine Mouse Antibody Isotyping **Sample Buffer** (Order No. MQSBXX, dep. on type):

Product Name	Order No.	Content	Preparation	Storage	Shelf life
Anti-mouse IgG1 Sample Buffer	MQSBG1	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse IgG2a Sample Buffer	MQSBG2A	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse IgG2b Sample Buffer	MQSBG2B	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse IgG2c Sample Buffer	MQSBG2C	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse IgG3 Sample Buffer	MQSBG3	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse IgM Sample Buffer	MQSBM	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse IgA Sample Buffer	MQSBA	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse kappa light chain Sample Buffer	MQSBKLC	0,5 mL	ready to use	2°C to 8 °C	until expiry date
Anti-mouse lambda light chain Sample Buffer	MQSBLLC	0,5 mL	ready to use	2°C to 8 °C	until expiry date

Additionally Required Materials

Important:

In addition to the Universal Module (essential, Order No. MQUM 1) and the individually selected Sample Buffer(s) (Order No. MQSBXX, dep. on type), the following items are required:

- 96-well microtiter plate
Pipets (2-20 µL, 20-200 µL) and tips

Intended Use

The Milenia QuickLine Mouse Antibody Isotyping System, consisting of the Universal Module (Order No. MQUM 1) and Sample Buffer(s) (Order No. MQSBXX, depending on type), is designed for fast and easy identification of mouse antibody isotypes, subtypes and light chains.

For research use only. Not for diagnostic purposes.

Principle of the Test

Knowing the antibody subclass is important when mouse monoclonal antibodies are produced. Hereby the optimal purification method can be chosen for the specific isotype which results in the highest yield of the antibody. Furthermore if there is no antigen available after recloning, positive clones can easily be detected with this test and monoclonality can be verified.

Milenia QuickLine Mouse Antibody Isotyping is a lateral flow immunoassay designed for the determination of mouse monoclonal antibody isotypes, subtypes and light chains within 5 minutes.

The immunoassay can be used for cell culture supernatants and/or purified antibodies.

First the isotype specific Sample Buffer (e.g. anti-mouse IgG1) is pipetted into a well of a 96-well microtiter plate. Subsequently the test sample of unknown antibody subclass is added. If a complex is formed between the Sample Buffer and the test sample, binding to an anti-mouse antibody on the universal test strip will result in a red line formation.

Test Procedure

Important notes:

All reagents must be equilibrated to room temperature (18-28°) before use.

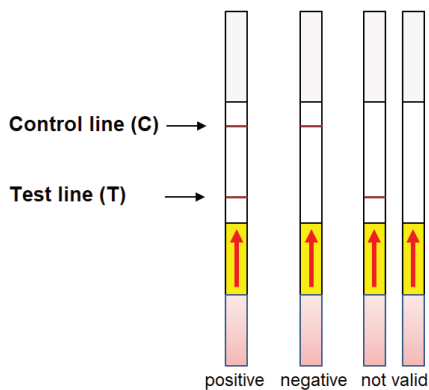
If the antibody concentration is too high, it can lead to a false negative result (high dose hook effect). This may occur when highly concentrated purified antibodies are applied in the test. In this case, samples should be prediluted with 1 X PBS to enter in the working range of the assay (1 – 20 µg/mL).

Test Setup:

1. Pipet **45 µL** of the appropriate **isotype-specific Sample Buffer (MQSBXX)** into a well of a 96-well microtiter plate
2. Add **10 µL** of the **antibody test solution** (about 1-20 µg/mL) and mix well
3. Place one **universal test strip** of the Milenia QuickLine Mouse Antibody Isotyping Universal Module (MQUM 1) into every well
4. Incubate for **5 minutes** at room temperature
5. **Read the results by visual inspection** immediately after incubation time

Note! Test lines developed after this time are without relevance

Interpretation of Results

Positive:

Test line (T) and control line (C) are visible

Specific isotype has been detected (valid pos result)

Note: faint red test line should be reported as positive

Negative:

Only control line (C) is visible

No isotype can be detected (valid neg result)

Not valid:

Only test line (T) or no line visible

Tests are interpreted visually. Examples of test results are shown in the figure above.

The **control line (C)** should be positive in all cases, indicating a successful test.
The **test line (T)** is associated with the presence or absence of the target isotype.

Troubleshooting

Problem	Possible cause(s)	Recommendation
No visible control line	a) damaged test strip	a) repeat test with new test strip
False negative result	a) concentration of purified antibody is too high b) concentration of antibody in cell culture supernatant is too low	a) dilute antibody in 1 X PBS b) use higher amount of antibody
Several test lines occur	a) incubation time was exceeded b) concentration of purified antibody is too high c) more than one antibody present in the sample	a) do not incubate for more than 5 minutes b) dilute antibody in 1 X PBS c) reclone cell line

If no line is visible on the test strip, please contact the Milenia Biotec GmbH customer service.

Assay Sensitivity

Sensitivity varies based on different conditions (purified antibody/cell culture supernatant). Antibody amounts between 1 and 20 µg/mL are detectable.

Documentation of Results

For the purpose of documentaion, it is possible to cut der lower part of the test strip (above the yellow area) after incubation time and, if desired, make photos of the strips.

Note: During drying process, the color of the result lines intensify.

Quick Reference Guide

1. 45 µL isotype specific Sample Buffer
2. 10 µl antibody sample, mix
3. place 1 test strip into well
4. 5 min incubation and read results immediately afterwards

For any questions and suggestions, please contact us:

Telefon: 0 641 – 94 88 83 – 0
Fax: 0 641 – 94 88 83 – 80
E-mail: info@milenia-biotec.de
Web: www.milenia-biotec.de

Mouse Antibody Isotyping Anleitung

Instruction for the Identification of Mouse Monoclonal
Antibody Iso-, Subtypes and Light Chains

Required components:

Mouse Ab Isotyping Universal Module (Order No. MQUM 1)

Mouse Ab Isotyping Sample Buffers (Order No. MQSBXX)

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Ist auf dem Teststreifen keine Linie sichtbar, wenden Sie sich an den Kundendienst der Milenia Biotec GmbH.
Milenia Biotec GmbH · Versailler Str. 1 · 35394 Gießen, Germany